

## Flow Cytometry Wash buffers

<b>PBS/BSA/Azide (PBA)</b>
1 litre sterile PBS 2 mL 20% sodium azide (final 0.04%) 10 g BSA (final 1%)

<b>PBS/Azide</b>
1 litre sterile PBS 2 mL 20% sodium azide (final 0.04%)

<b>PBS/AB/Azide (PBA)</b>
1 litre sterile PBS 2 mL 20% sodium azide (final 0.04%) 10 mL heat-inactivated AB serum (final 2%)